D

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thermocult

Operating Instructions

Small-Size Incubator with permanently set incubation temperature


- permanently set incubation temperature of 35°C +/-2°C
- economical space configuration with horizontal or vertical mounting or placement. Suitable for 20 dip-slide cultures, 8 Petri- dishes or 6 blood culture bottles.
- 2 separate control lights for power and heating system
- built in thermometer
- Shelf for tray cultures and staining racks/dishes (optional)
- All-around heating elements for uniform heat distribution
- Clear door-visual control, rapid check
- Radio interference suppression Degree N
- Casing of high quality poly-material
- Common cleaners and disinfectants can be used

Table of Contents

Page

| | |
|-----------------------------------|----|
| Control elements | 9 |
| Placement of the unit | 9 |
| Start-up | 10 |
| Working temperature | 10 |
| Incubation temperature adjustment | 10 |
| Hints for the user | 10 |
| Cleaning the incubator | 10 |
| Troubleshooting | 10 |
| Technical specs. | 11 |
| Statement of conformity | 12 |

Control Elements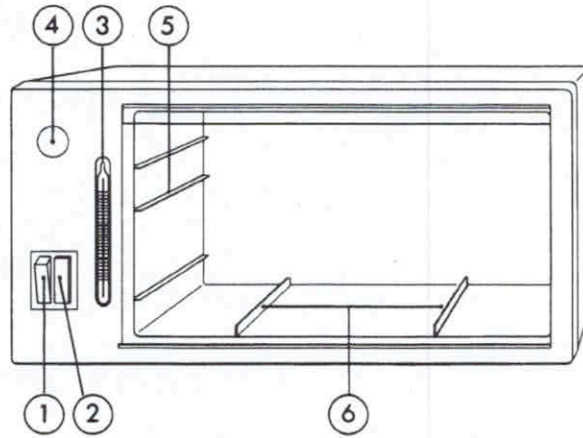

- 1 Master switch with Service control light (down: ON)
- 2 Heating System control light
- 3 Thermometer
- 4 Cover for the temperature adjustment screw
- 5 Support ledge for shelves
- 6 Bottom Strips for slide cultures

Accessories (not shown): shelves

1. Placement of the unit

The Small-size incubator thermocult can be used in a horizontal (Fig.1) or a vertical (Fig. 2) Position.

When setting the incubator upright, make sure to attach the foam rubber feet. The incubator **should not sit directly on the pad**

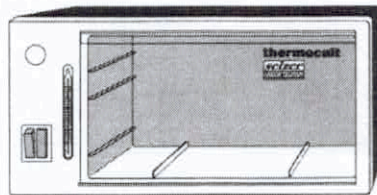
Fig. 1

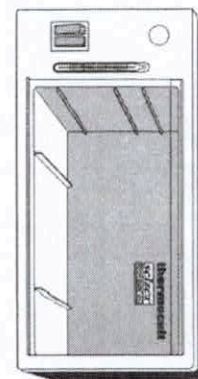
Fig. 2

2. Start-Up

- a) Check to ensure that the voltage indicated on the rating plate and the Power supply voltage match.
- b) Plug the line plug into an electrical outlet (230V ~)
- c) Turn on the master switch (1): The red Service control light (1) and the yellow heating System control light (2) are on.
- d) The constant chamber temperature is reached after approx. **60 minutes**. The yellow heating System control light (2) will turn on and off at irregular intervals.

3. Working Temperature

The built-in temperature controller maintains a constant Incubation temperature of 35°C with a variation of $\pm 2^\circ\text{C}$ at various spots of the chamber (ambient temperature: 20°C - 25°C).

4. Incubation Temperature Adjustment

An adjustment is required if, after prolonged use of the Instrument, the permissible fluctuation span is no longer maintained. In order to proceed, remove the cover(4). If you turn the screw behind it to the right, you can adjust a higher incubation temperature. You can lower it by turning the screw to the left.

A one-quarter rotation corresponds to approx. 15°C.

5. Hints for the user

To close properly, the door must be pressed shut until impact. The thermocult unit must be installed in a Position where the door is not exposed to direct sun light.

The shelf should be used for all vessels or plates **so that they do not sit directly on the floor.**

It is recommended to switch off the device at longer disuse.

6. Cleaning the Incubator

Both the inner and the outer casing as well as the door of the incubator consist of high quality poly-materials.

The surfaces will tolerate all commonly used cleaners and disinfectants.

Solvents (e.g. methanol, acetone, and the like) should not be used to clean the unit.

7. Trouble-shooting

Have the unit returned to the manufacturer by your dealer. If the Problem falls within the warranty, make sure to include the warranty certificate and add a short description of the fault.

Technical Specs.

| | |
|-------------------------------|---|
| Type | thermocult |
| Temperature constancy | $\pm 0.3^{\circ}\text{C}$ |
| Incubation temperature | $35^{\circ}\text{C} \pm 2^{\circ}\text{C}$ |
| Rated voltage | 230V ~ |
| Current and frequency | A.C. ~50/60 Hz |
| Rated current | 0.14 A at 230 V |
| Rated capacity | 30W |
| Class of protection | II (protective insulation) Not "ex" proof |
| Radio interference suppressed | degree N (VDE0875) |
| Technical safety compliance |  |
| Capacity | 20 dip-slide cultures, or 8 Petri dishes or 6 blood culture bottles |
| Dimensions: | |
| Casing | 30 x 17 x 15cm (l x d x h) |
| Work chamber | 21 x 14.5 x 11cm (l x d x h) |
| Weight | approx.1.3kg |
| Accessories | shelves |

**KONFORMITÄTSERKLÄRUNG / DECLARATION OF CONFORMITY /
DECLARATION DE CONFORMITE/ DICHIARAZIONE DI CONFORMITA**

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Meets all the provisions of the directive **2014/35/EU; 2014/30/EU; 2011/65/EU** what apply to it.

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